



VICON NANTAGE

QUICK START GUIDE

Contents

Check what's in the box 3

Set up your Vicon Vantage system 5

Configure your Vicon Vantage system 9

Vicon Vantage troubleshooting 12

© Copyright 2015–2020 Vicon Motion Systems Limited. All rights reserved.

Vicon Motion Systems Limited reserves the right to make changes to information in this document without notice. Companies, names, and data used in examples are fictitious unless otherwise noted. No part of this publication may be reproduced, stored in a retrieval system, or transmitted in any form or by any means, electronic or mechanical, by photocopying or recording, or otherwise without the prior written permission of Vicon Motion Systems Ltd.

Vicon® is a registered trademark of Oxford Metrics plc. Vicon Bonita™, Vicon Lock™, Vicon Lock+™, Vicon Lock Lab™, Vicon Lock Studio™, Vicon Nexus™, Vicon MX™, Vicon Shōgun™, Vicon Tracker™, Vicon Vantage™, Vicon Vero™, and Vicon Vue™ are trademarks of Oxford Metrics plc.

VESA® is a registered trademark owned by VESA (www.vesa.org/about-vesa/). Other product and company names herein may be the trademarks of their respective owners. For full and up-to-date copyright and trademark acknowledgements, visit <https://www.vicon.com/vicon/copyright-information>.

Vicon Motion Systems is an Oxford Metrics plc company.

Email: support@vicon.com Web: <http://www.vicon.com>

<https://accessgudid.nlm.nih.gov/>

B548VICONVANTAGESYSTEM0



Check what's in the box

Before setting up your Vicon Vantage system, check that you have all the necessary components, which are described in this section.

- ✔ For a list of components supplied with your specific Vicon system, see your shipment Bill of Materials.

Hardware

Vicon Vantage cameras (V16, V8, V5)

Vicon optical cameras with custom-engineered sensors and lenses, high-power strobes, on-board processing, Gigabit Ethernet connectivity and intuitive design. Available in high-performance resolutions up to 16 megapixels at 120 fps full-frame capture speeds.



PoE+ switch

The supplied PoE+ switch provides the power source and connectivity for your Vicon Vantage system.

Vantage cameras require a PoE+ switch.

Vicon Lock units (via a PoE switch) and Vicon Vero and Vue cameras can also connect to a PoE+ switch.



Vicon Lock unit

Vicon's optional connectivity device for the Vantage system. The Vicon Lock unit (such as the Lock+, Lock Studio, or Lock Lab) synchronizes with your Vantage, Vero and Vue cameras. It also provides interfacing to third-party analog devices, timecode, genlock and sync out functionality.



Vantage host PC

A high-performance Vicon-supplied PC, enabling communications between the PoE+ switch and the installed Vicon application software. This is where you capture, view, and process the data from the Vantage system and any supported third-party devices integrated into it (Advanced spec PC illustrated).



Vantage cables

You can connect up a simple Vantage system consisting of 1–12 Vantage cameras and a PoE+ switch, to the Vicon Vantage host PC, using the supplied RJ-45 cables. Camera cables include a ferrite at the camera end only. If you are using one or more Vicon Lock+ units, or MX T-Series cameras with a Gigaset, additional cables are supplied.



Vantage calibration device

Vicon's Active Wand calibration device makes it quick and easy to calibrate the Vantage cameras as well as video cameras, and your capture volume.



Software and accessories

Vantage accessories

Vicon motion capture consumables include marker fixing tape, arm and head bands, anthropometric measuring tools, Micropore and gaffer tape, Velcro, unique Vicon retro-reflective markers in a variety of sizes, Lycra® suits, and more. The consumables that are supplied depend on which items you've ordered with your Vicon system.



Product installation download

Your Vicon Vantage installation files, available on the Vicon website, include your licensed Vicon application software and VAULT licensing software.



www.vicon.com/downloads

Set up your Vicon Vantage system


Define the capture volume

- 1 Place markers around the edges of the volume to indicate the intended capture volume.

You can often place the markers on the floor, as shown below. If the volume to be captured is higher up, place the markers accordingly. This makes it easier to visualize the capture space.

- 2 Roughly position the cameras around the volume to provide full coverage across the volume.

You can find more information in [Position the cameras \(page 8\)](#). Positions will depend on the volume size and camera lenses.

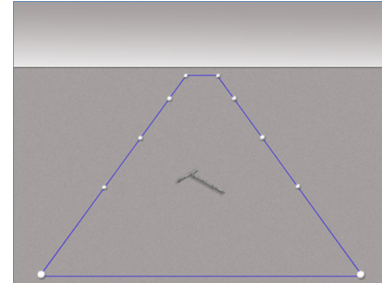

- ✓ Remember to include height as well as width and depth in your volume calculations!

Install and connect the cameras

- 1 Install each camera on a tripod, on the wall, or on a ceiling-mounted rig.

You can order mounting brackets and camera mounts from Vicon.


- 2 Ensure the Power over Ethernet (PoE+) switch is connected to an appropriate power source.
- 3 For each camera (up to 12 per PoE+ unit):

- Insert the ferrite end of one of the supplied network cables (RJ-45) into the camera.
- Connect the other end of the cable to the PoE+ switch.



Although there are 26 sockets in total, always remember:

- One socket is needed to connect to the Vicon Vantage host PC.
- The 24 PoE ports can support up to 12 Vantage cameras and any of the other Vicon components in your system, such as a Vicon Lock unit (connected via an additional PoE+ switch), Vicon Vue video cameras, etc. If required, the uplink ports or any remaining PoE ports can be connected to the PC.

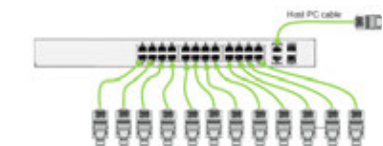


To add more cameras, you can connect further PoE+ switches to the system (see Vicon system configuration and connection examples, or the Vicon Vantage Reference Guide).



- 4 Turn on the PoE+ switch.


Connect and configure the PC

- 1 With the supplied network cable (RJ-45), connect the PoE+ switch to the PC.
- 2 On the PC open **Windows Network Connections**, right-click the NIC to which the system is to be connected and then click **Properties**. (If you are connecting more than one card, right-click the first NIC.)
- 3 In the **Properties** dialog box, on the **Networking** tab, clear all the unwanted items, leaving only **Internet Protocol Version 4 (TCP/IPv4)** selected.

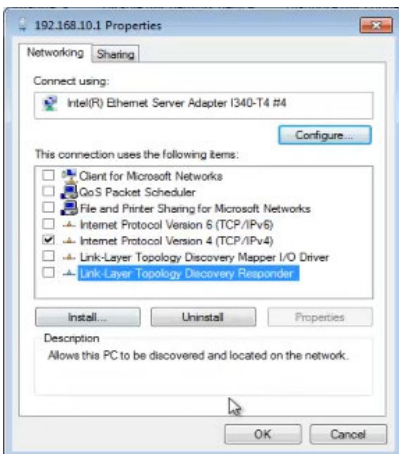

✓ For detailed instructions on configuring your PC network card, see [Configuring network card settings](#)¹ or *Configuring and connecting the network ports* in the Vicon PDF, *PC Setup for Vicon systems*.

- 4 Click on this option to highlight it and then click **Properties**.
- 5 In the **Internet Protocol Properties** dialog box, on the **General** tab enter the following information:
 - **IP address** box, enter 192 . 168 . 10 . 1.
 - **Subnet mask** box, enter 255 . 255 . 255 . 0.
 - **Default gateway** box, leave the box blank.

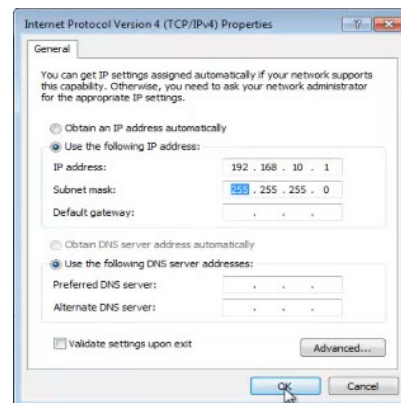

- 6 If you are using more than one NIC, enter the **IP address** 192.168.10.1 for the first NIC to which the system is connected, as described above, see the following *Link aggregation for multiple NICs* box for the remaining addresses.
- 7 Click **OK** to close the dialog box, and **OK** to close the Properties dialog box and save the settings.

¹ <https://docs.vicon.com/display/Connect/Configuring+network+card+settings>

✔ Link aggregation for multiple NICs

In very large systems, to enable you to use more cameras or data, you can use additional NICs in your PC to split camera data. To do this:

1. Set the additional IP addresses of your NICs to different addresses in the available range 192.168.10.2–9.
2. In your Vicon application software, in the Properties pane for each camera, set the **Destination IP Address** to the IP address of the NIC to which to send data.

Install the Vicon application software

1 Download the required software installer from the Downloads page on the Vicon web site www.vicon.com/downloads².

2 Double-click on the installer (setup.exe), read and accept the User License Agreement, and follow the instructions on each page.

Your Vicon software automatically installs into its own folder.

The Vicon software installer enables optional installation of the License Server software for server-based licensing, and the dongle drivers for licenses that are locked to dongles.

3 If you are prompted to restart your PC, do so as soon as possible.

License the Vicon application software

Vicon application software uses Safenet licensing and a built-in licensing utility (the Vicon Automated Unified Licensing Tool, VAULT).

You specify the license server in a dialog box that appears if no license is found or via the **Help > Vicon Product Licensing** menu option. In Windows 10, the server name is stored in the following file:

```
<InstallationDrive>:\Users\Public\Documents\Vicon\Licensing\<ProductName>LicInfo.xml
```

If you need to view information about licensing while you are using your Vicon software, on the **Help** menu, click **Vicon Product Licensing**.

For detailed instructions on how to use VAULT, see your Vicon software documentation (PDF or online help).

² <http://www.vicon.com/downloads>

Start the software




1 On your Windows desktop, double-click the shortcut for the required Vicon software.



2 If this is the first time you have used your Vicon application software on this PC, a dialog box may be displayed, enabling you to specify the license server.

3 In the System Resources pane, make sure all of the cameras in your system are listed.


4 Ensure that at the top left of the screen the system is shown as live and connected.

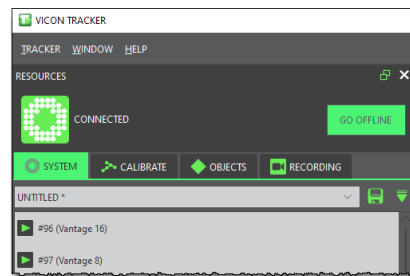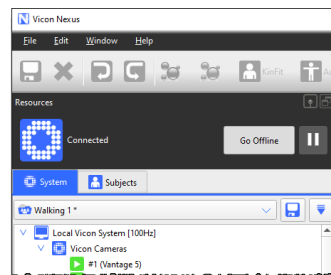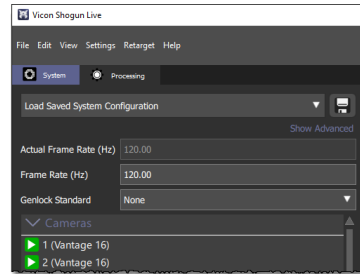

Position the cameras

- 1 In the System Resources pane, select a camera.
- 2 Change the View pane to Camera view.
- 3 Aim the camera so that it points where required, usually the center of the capture volume.
- 4 Repeat the process for each camera.

The diagram (right) shows typical camera coverage of the capture volume.

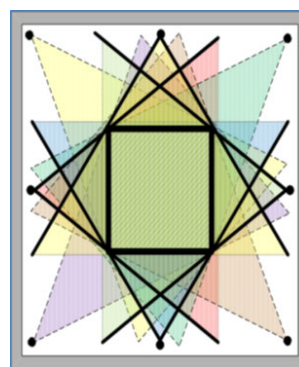

Configure your Vicon Vantage system

Adjust camera settings

- 1 Ensure you are in Live mode and that a **Camera** view is displayed.
- 2 In the **System Resources** pane, select one camera.
- 3 Notice that the **Properties** for this camera are listed at the bottom of the pane.
- 4 Have someone wave the calibration wand in the capture volume and ensure that you can see markers moving in the view pane.
- 5 Note the **Grayscale Mode** property. The default setting is Auto, which tells the cameras to send to the Vicon software all the centroid data for grayscale blobs it can circle fit. If any grayscale blobs cannot be circle fitted, the grayscale data is sent to the Vicon software so it can try to circle fit the blobs.
- 6 Look at the **Strobe Intensity** setting. With a value of 1, the cameras send as much light as possible into the volume. Zoom in to a marker by right-clicking and dragging the mouse. Notice how the pixels change as you lower the light (to view pictures from a grayscale blob, **Grayscale Mode** must be set to All).
- 7 Notice the **Threshold** property. It tells the camera how bright a pixel must be to be accepted by the camera. Zoom in to a marker and notice how the pixels change as you adjust the property.
- 8 Adjust the **Strobe** and **Threshold** settings so that you see only the 2D circular-shaped markers in your volume. Adjust the **Strobe Intensity** setting only as low and the **Threshold** only as high as needed.

Set camera focus and aperture

- 1 Place three markers in the middle and on the far side of the capture volume.
- 2 Leaving the focus on infinity (∞), close the aperture to as high an f-stop value as possible, so that you can still see the markers at the furthest distance required.

✓ In general, it's better to run with a low Threshold setting and more closed aperture, than a high Threshold and open aperture. (See **Threshold** in [Adjust camera settings \(page 9\)](#).)



The illustration (right) shows the focus and aperture rings on a Vicon 18 mm lens.

3 For each camera, turn the aperture ring (the ring with the f-stops) until the aperture is fully open. Later you will turn the ring back, increasing the f-stop number and closing the aperture.

4 In the **System Resources** pane, select all the cameras by Shift-clicking. In the **Properties** section at the bottom of the window, change **Grayscale Mode** to All.

5 In the **System Resources** pane, select a camera. For each marker:

1. In the view pane, zoom in to the center of the marker.


⚠ Vantage cameras are supplied pre-focused for typical motion capture use, so in most situations, no focusing is necessary. The following step (b) is applicable only if you need to re-focus the camera, otherwise, go straight to step c.

2. If required, turn the camera focus ring (shown in the previous illustration) until the 2D image of the marker in the view pane is clear and sharp. As you rotate the focus ring, note that the marker images get smaller and then begin to get larger again. Good focus is when the marker is at its smallest size.
3. Observe that, with the aperture fully open, the markers appear fully saturated, with very little grayscale gradation and some merging, as in the following image.



Adjust the aperture ring to change the amount of light that hits the image sensor. Turn the aperture ring to close the aperture and obtain good, clear images of markers even when they are close together.

If you turn the aperture ring too far, until the aperture is almost closed, the markers appear very small, lack any bright grayscale pixels and the circle-fitting is poor, as shown below.

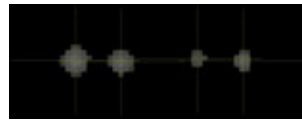

Adjust the aperture ring so that the markers are not too small, and the center is just off-white (not fully saturated) when viewed from the middle of the volume.



6 Repeat Step 5 for each camera.

7 When you are happy with the focus and aperture, to prevent the rings from moving if the camera is knocked, tighten the locking screws.

⚠ To avoid damaging the rings, do not over-tighten the locking screws.

8 Select all of the cameras and change the **Grayscale Mode** property back to Auto.

Mask unwanted reflections

- 1 Ensure that your camera settings are as recommended in [Adjust camera settings \(page 9\)](#).
- 2 From the capture volume, remove any unnecessary objects, such as the calibration device, markers, or other objects likely to cause reflections.
- 3 Make sure **Camera** view is displayed. In the **System Resources** pane, select all cameras.

- 4 In your Vicon software, click **Start**:
 - Nexus: **System Preparation Tools** pane > **Mask Cameras** section
 - Tracker: **Calibrate** tab > **Create Camera Masks** section
 - Shogun Live: **Camera Calibration Tools** pane > **Wand** section
- 5 After a few seconds, when all reflections appear to have been covered, click **Stop**.
- 6 If you still see unwanted reflections, you can start and stop camera masking again until you are satisfied. Alternatively, use the **Manual Camera Masking** tools.

Calibrate the cameras


You calibrate the cameras to enable the cameras to capture subject movements throughout the volume.

You can also set the volume origin and orientation of axes in relation to the capture volume and to one another.

To calibrate your Vicon Vantage cameras, use the supplied Vicon calibration device.

- 1 Remove from the capture volume any unnecessary objects, such as calibration objects and markers.
- 2 In your Vicon application software, ensure the following are selected:
 - All the cameras
 - The calibration device you will be using
 - An appropriate number of **Refinement Frames**
 - **Auto Stop**
- 3 In the **Calibrate Cameras** section, click **Start**.

- 4 In the volume, wave the wand to cover the whole target volume area. Ensure that the tracks are evenly distributed throughout the volume. Check the distribution of the tracks in the view pane.

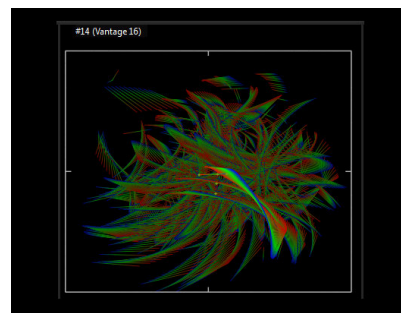

When all cameras have captured the number of frames set by the Refinement Frames option, the calibration stops.

Additional information about setup, calibration and object tracking can be found in the online help of your Vicon software.

Vicon Vantage troubleshooting

If something is not quite right, try these quick tips to get things going again.

What's in the box

I am missing some equipment
or

Some equipment appears damaged from shipment.

- Contact Vicon Support support@vicon.com³

Connecting

My hardware isn't connecting to the computer.

- Confirm that the IP address settings are correct following the instructions in [Connect and configure the PC \(page 6\)](#).
- Check the network card settings, including Jumbo Packet (for details, see *Configuring network card settings* or *Configuring and connecting the network ports* in the PDF *PC setup for Vicon systems*).
- Disable any firewall on the network card.
- Confirm that the Ethernet cable is firmly seated.
- Ensure you are an administrator on the computer.

Camera hardware

I don't see an image from the camera.

- Confirm that you've removed the lens cap.
- Ensure that the aperture is not fully closed.
- Confirm that the **Strobe Intensity** is high (best on full at 1) and that the **Threshold** is not too high (a good default is 0.2).

When I tap a calibrated camera to select it, the camera indicates that it has been bumped.

- To change the sensitivity of the camera to taps, in the camera's **Properties** pane, **Settings** section (Nexus and Tracker) or **General** section (Shogun Live), choose a lower **Bump Sensitivity** setting.

Calibrating cameras

After calibrating, image errors are very high.

- Ensure that the camera aim and focus are correct, the camera properties are set appropriately, and that you covered the whole volume during your wand wave, then repeat the calibration.
- Ensure **Active Wand** is selected in the **Calibrate Cameras** settings (Nexus and Tracker) or **Camera Calibration** tab (Shogun Live).

If you need more information than that supplied in the documentation or on the [Vicon Support web pages](#)⁴, please contact Vicon:

Denver, CO
Vicon Denver
7388 S. Revere Parkway
Suite 901, Centennial
CO 80112, USA
T: 303.799.8686
F: 303.799.8690
E: support@vicon.com

Los Angeles, CA
Vicon LA
3750 S. Robertson Boulevard
Suite 100, Culver City, Los Angeles
CA 90232, USA
T: 310.437.4499
E: support@vicon.com

Oxford, UK
Vicon Oxford
Unit 6, Oxford Industrial Park
Mead Rd, Yarnton, Oxford
OX5 1QU, United Kingdom
T: +44.1865.261800
E: support@vicon.com

³ <mailto:support@vicon.com>

⁴ <https://www.vicon.com/support>