

# Capture.U


## QUICK START GUIDE



What's in the box?  
Blue Trident sensor(s)  
Sensor strap(s)  
IMU adapter(s)  
Micro USB cable(s)



### 1. Get to know the sensor

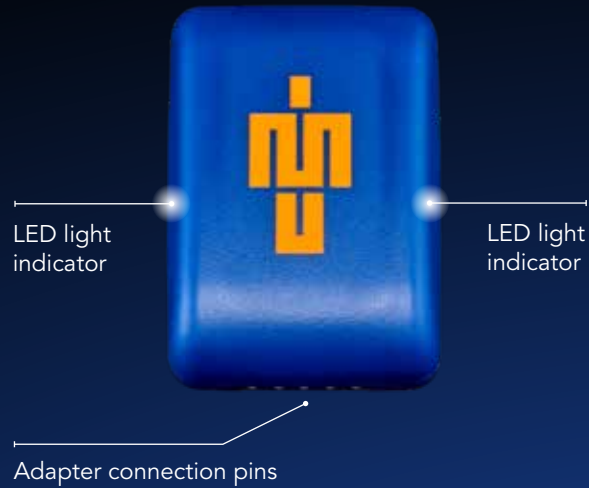

### 2. Download the Vicon Capture.U app from the App Store to your iOS device




### 3. Select a capture mode

- **To Sensor:** Capture data to sensor on-board memory (more than 6 sensors).
- **To Device:** Capture data to iOS device (up to 6 sensors).
- **Real-Time Insight:** Stream real-time data to iOS device with video overlay (2 sensors).
- **Activity Widgets:** Select a focused workflow for specific activities.




Tap to explore the Real-Time Insight mode.


### 4. Tap Select Sensors (bottom left)

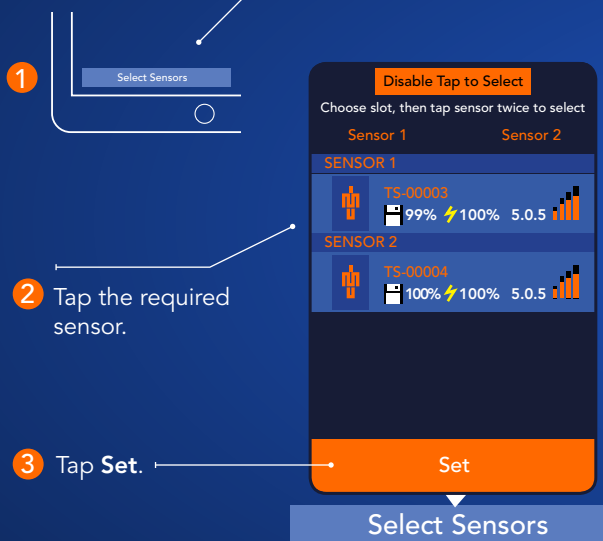


### 5. Select the data to stream and change the axis graph

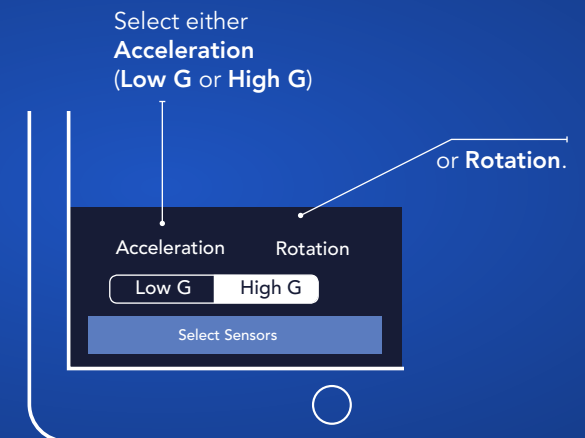

## 6. Tap the Video icon (top right)

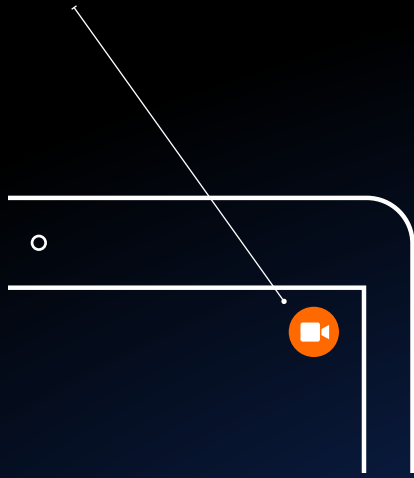


## 7. Set the benchmarks (bottom right)

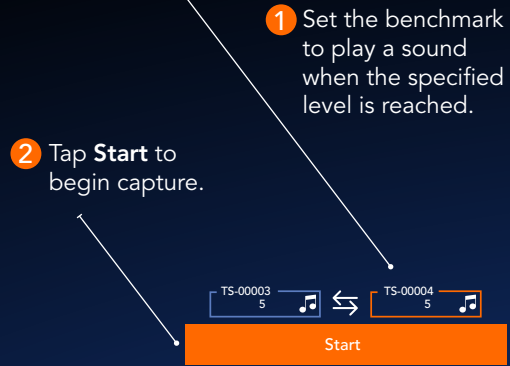


## 8. Tap the Statistics icon (bottom right)


When finished, tap **Stop**.

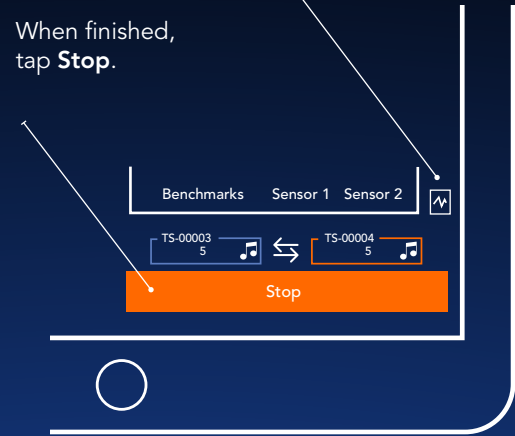

## 9. Select an export option:

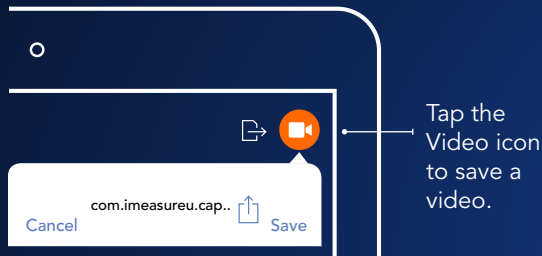


## 10. Tap Export to PDF to display a PDF Preview

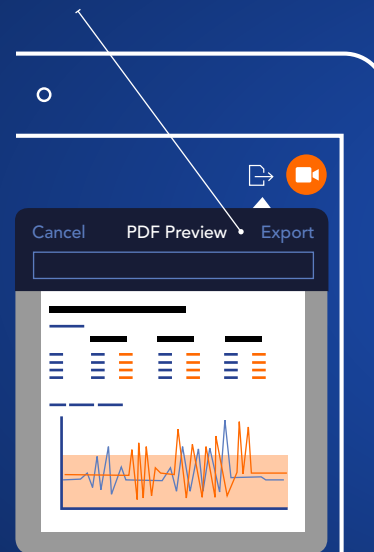

## 11. Download Capture.U Desktop [www.vicon.com/downloads/software](http://www.vicon.com/downloads/software)

Analyze your data and access walk-throughs and documentation.

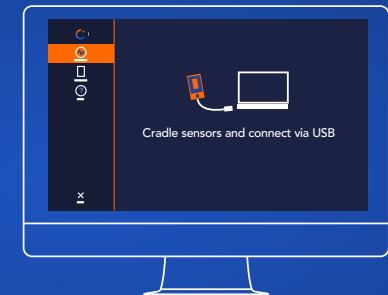

For regulatory information and links to other language versions of this guide, go to [docs.vicon.com/display/IMU/](http://docs.vicon.com/display/IMU/)

Copyright © 2019 Vicon Motion Systems Ltd. All rights reserved. Vicon® is a registered trademark of Oxford Metrics plc. IMeasureU is a trademark of IMeasureU Ltd in New Zealand. Other product and company names herein may be the trademarks of their respective owners. Vicon Motion Systems is an Oxford Metrics plc company. IMeasureU is a wholly owned subsidiary of Vicon Motion Systems Ltd.

[www.vicon.com/bluetrident](http://www.vicon.com/bluetrident) [www.vicon.com/captureu](http://www.vicon.com/captureu)