

IMeasureU Box



IMU Sensors



Sensor Straps




Micro-USB cables

Start Here


VICON

IMeasureU

01 The Sensor

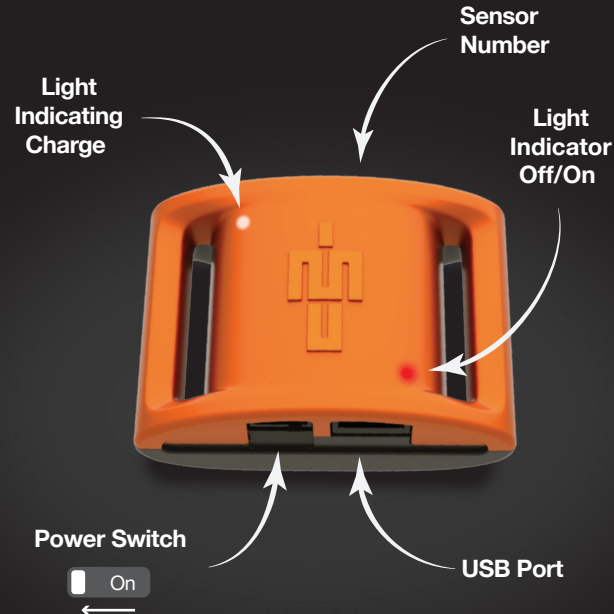


02 Download IMU Research App from the App Store with your iPhone or iPad

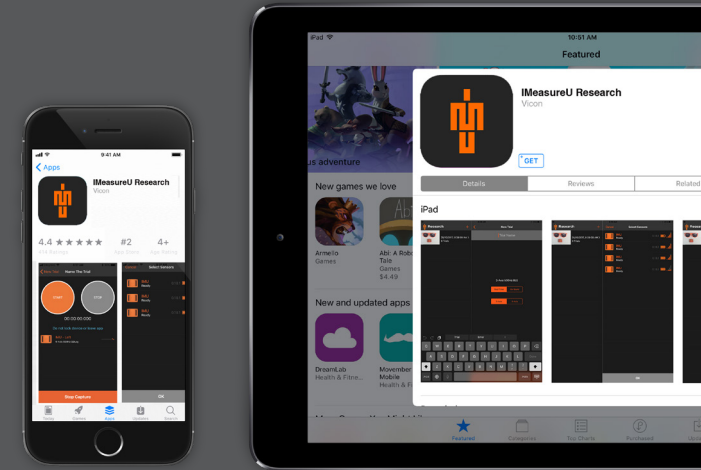


Unfold Me

03 Open the IMU Research App and create a new session

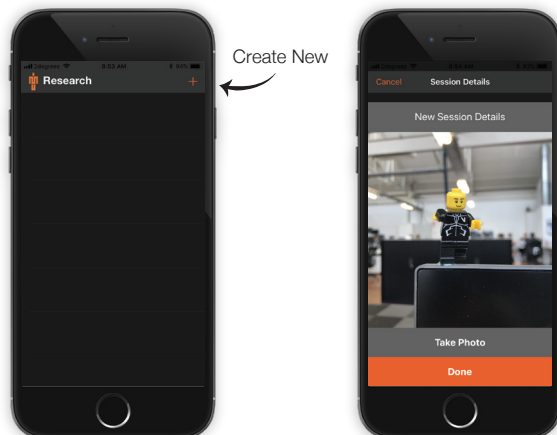


04 Create & record a new trial

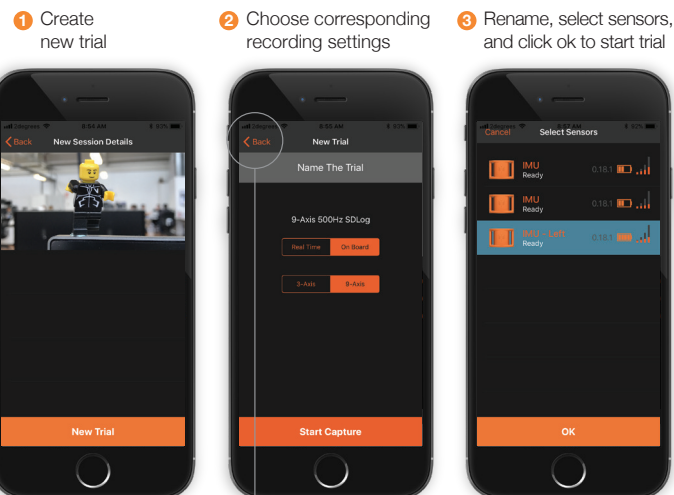

If you have multiple trials running, select 'back' to view all trials

05 Stop capture & export trial

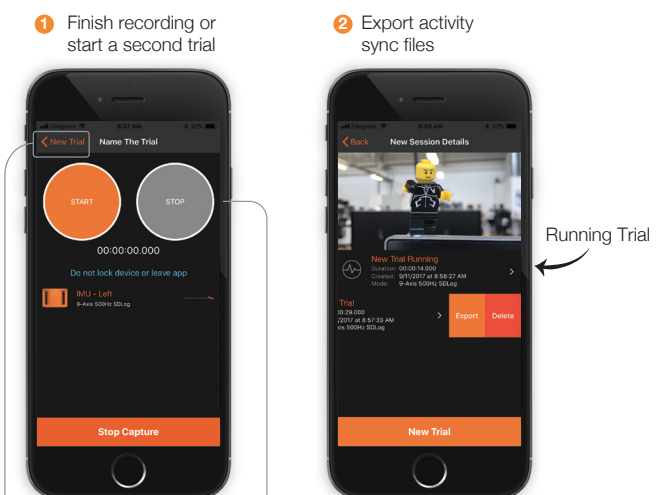

Start/stop buttons are to add user epochs into the sync file.

Flip Me

01 Make sure sensors are turned on before plugging in



02 Downloading preview data




03 Select and start download

1 Select desired sensor by serial number

2 Select exported sync file*

3 Select All /Add Regions of data

4 Download data

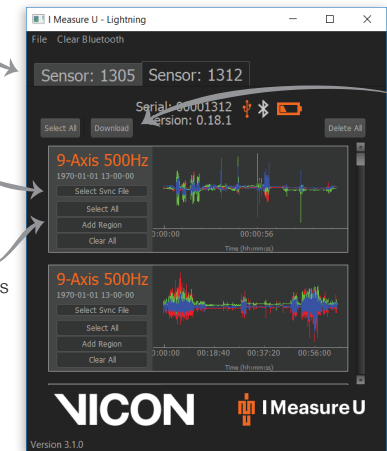

*Sync file text will turn blue once file is selected

04 Downloading and saving data


Saving and writing data*



*User will be notified of saving & verifying data when status bar is at 100%



05 Start new download or clear sensor

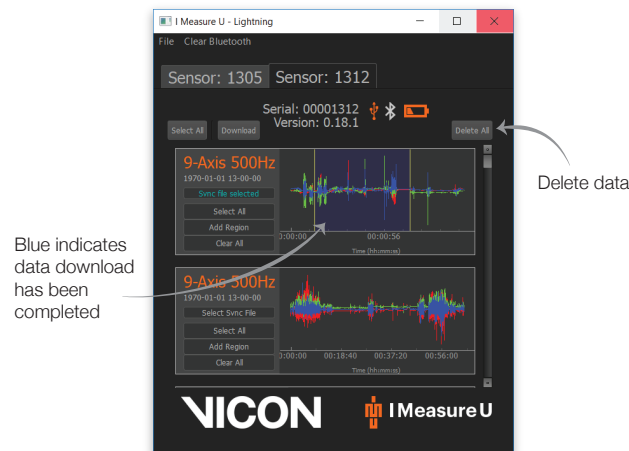

Back to step 3

For IMU regulatory information, see the enclosed document, Vicon IMeasureU Sensor V1 Safety and Regulatory Information.

© Copyright 2017 Vicon Motion Systems Limited. All rights reserved. Vicon® is a registered trademark of Oxford Metrics plc. IMeasureU is a trademark of IMeasureU Ltd in New Zealand. Other product and company names herein may be the trademarks of their respective owners. Vicon Motion Systems is an Oxford Metrics plc company.

www.vicon.com

www.IMeasureU.com

Thanks for reading us!

