

# Capture.U


## QUICK START GUIDE



What's in the box?  
Blue Trident sensor(s)  
Sensor strap(s)  
IMU adapter(s)  
Micro USB cable(s)



### 1. Get to know the sensor

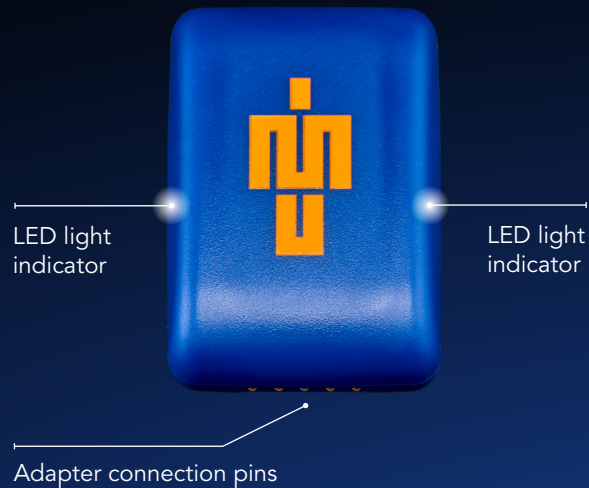

### 2. Download the Vicon Capture.U app from the App Store to your iOS Device



### 3. Select a mode:

Capture



TO SENSOR



TO DEVICE

Visualize



REAL-TIME INSIGHT



AR VISUALIZATION

Learn




EDUCATION




PRACTICE

Tap to explore the Real-Time Insight mode.


### 4. Tap Select Sensors (bottom left)

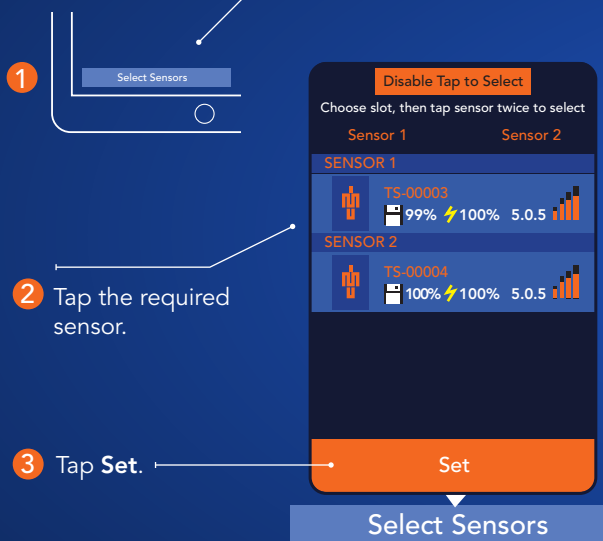

### 5. Select the data to stream and change the axis graph

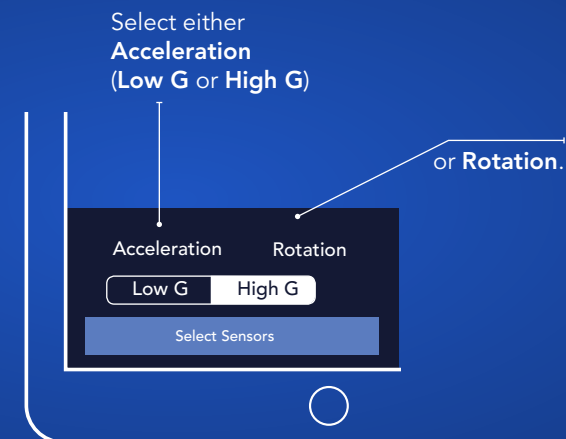


## 6. Tap the Video icon (top right)

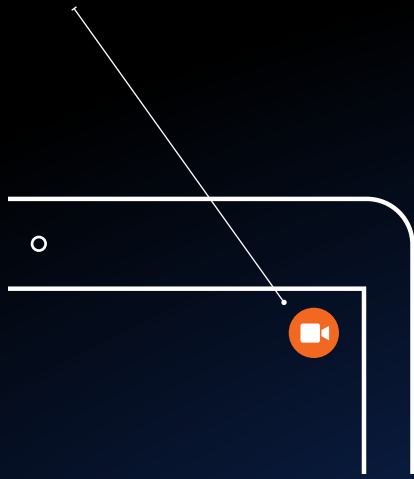


## 7. Set the benchmarks (bottom right)

- 1 Set the benchmark to play a sound when the specified level is reached.
- 2 Tap **Start** to begin capture.

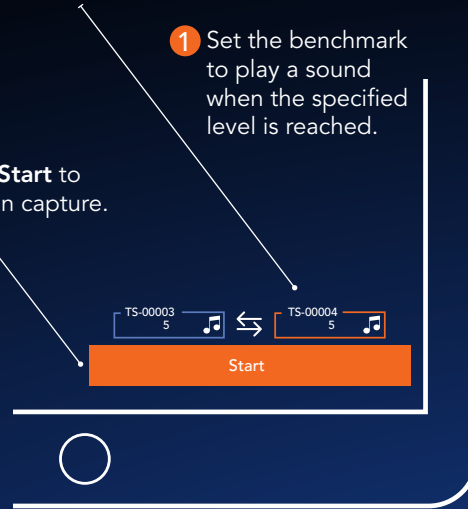


## 8. Tap the Statistics icon (bottom right)

When finished, tap **Stop**.

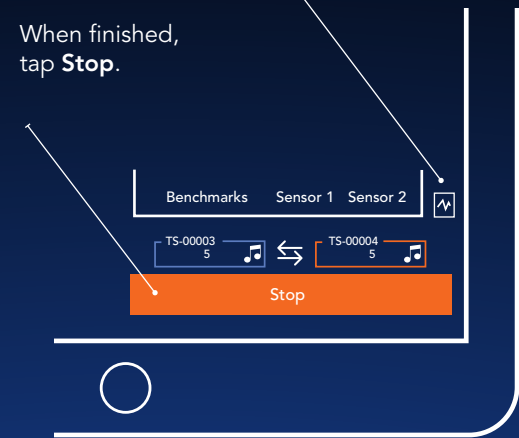

## 9. Select an export option:

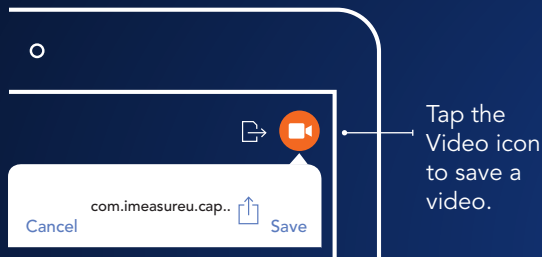


Tap the Video icon to save a video.




Tap the File icon to display the export to file options (PDF or CSV).

## 10. Tap Export to PDF to display a PDF Preview

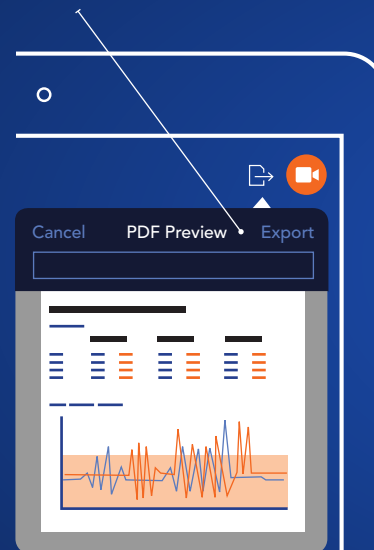

## 11. Download Capture.U Desktop [www.vicon.com/downloads/software](http://www.vicon.com/downloads/software)

Analyze your data and access walk-throughs and documentation.

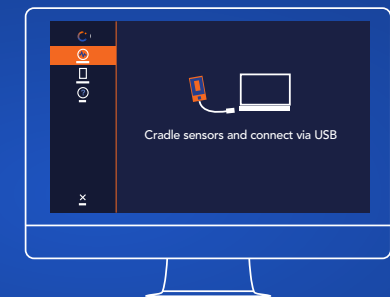

For regulatory information and links to other language versions of this guide, go to [docs.vicon.com/display/IMU/](http://docs.vicon.com/display/IMU/)

Copyright © 2019–2022 Vicon Motion Systems Ltd. All rights reserved. Vicon® is a registered trademark of Oxford Metrics plc. IMeasureU is a trademark of IMeasureU Ltd in New Zealand. Other product and company names herein may be the trademarks of their respective owners. Vicon Motion Systems is an Oxford Metrics plc company. IMeasureU is a wholly owned subsidiary of Vicon Motion Systems Ltd.

[www.vicon.com/bluetrident](http://www.vicon.com/bluetrident) [www.vicon.com/captureu](http://www.vicon.com/captureu)