

Release Notes Vicon Video Viewer 1.12.0

Vicon Motion Systems Ltd

Exported on 05-Jul-2023

Fixes

- Fixed bug performing distortion correction causing erroneous undistortion when principal point is not in image center.
- Added option (--centre-principal-point) to shift undistorted video such that the principal point is set to the image center.
- Fixed flipped video in ViconVideoConverter when processing H.264 video without changing format.
- ViconVideoConverter applies gamma correction and saturation when converting to H.264.
- UI remains responsive while large videos are opening in ViconVideoViewer.